

CSI ClorDiSys Solutions, Inc.

"The Chlorine Dioxide People"

Providing you with gaseous chlorine dioxide solutions for your decontamination and sterilization needs

Room Wall Plate

(Part #: CSI-ROOM008 with RH Probe)

(Part #: CSI-ROOM009 w/out RH Probe)

Description


The Room Wall Plate provides a method to connect ClorDiSys Solutions, Inc gas generators to a room for quick, easy connections. It is mounted through a studded wall to the room to be decontaminated.

The plate allows for quick valved tubing connections and quick connection of electrical connectors to measure and control the RH to either the Mindox or Cloridox systems



Assembly

1. Cut wall and drill holes as per Wall Cutout Template
2. Insert Gasket over protruded side of Wall Plate
3. Mount Gasket and Wall Plate into the wall, with the indented side with the connectors facing the outside of the Decon room.
4. Insert cement screws into through holes of Wall Plate and tighten.
5. Position Bezel Plate to the wall on the Decon room side of the assembly.
6. Caulk Bezel Plate in place.

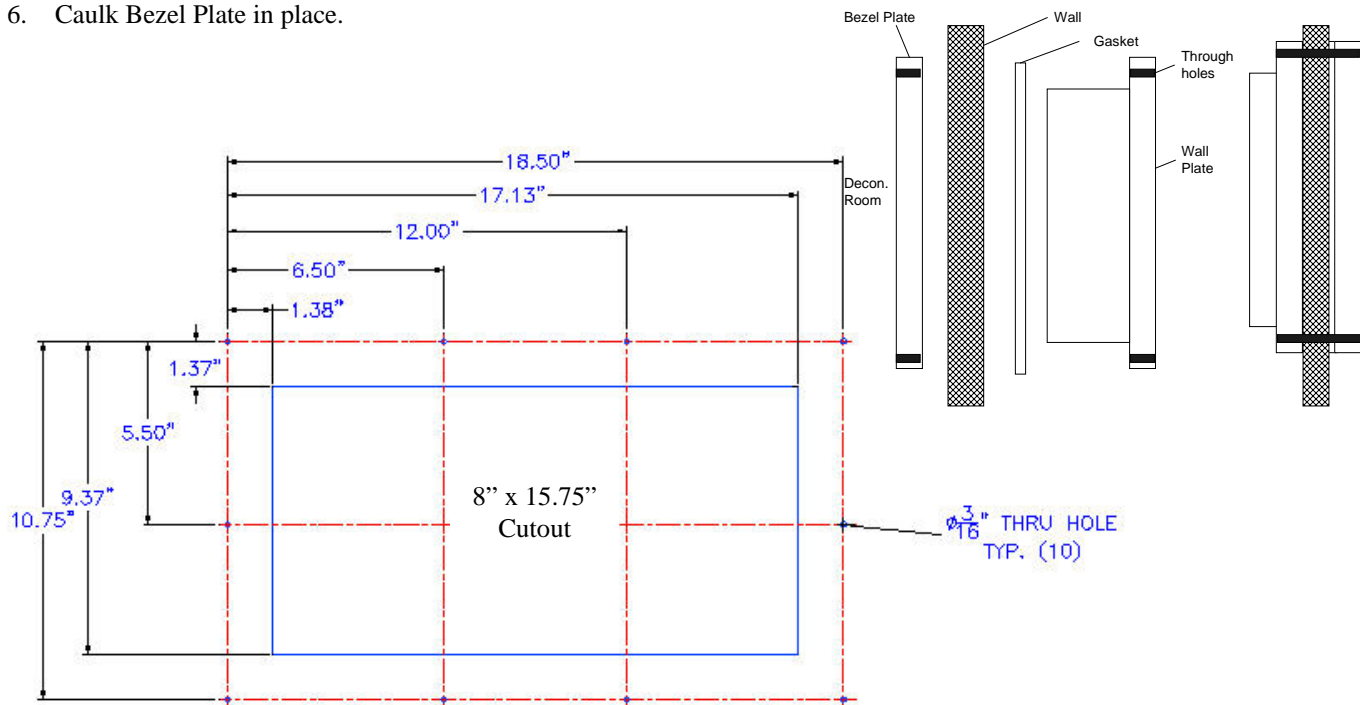

CSI ClorDiSys Solutions, Inc.

P.O. Box 549, Lebanon, NJ 08833-0549 Tel:(908) 236-4100 Fax:(908) 236-2222

www.clordisys.com

100-009 Rev B