



"The Chlorine Dioxide People"

Providing you with chlorine dioxide solutions for your decontamination needs

CD Check Strips

Description:

Get an easy, quick and cost-effective way to document your chlorine dioxide decontamination process. A CD Check Strip™ changes color when exposed to chlorine dioxide gas. The color change provides an indication that the strip was exposed to chlorine dioxide gas. This color change does not guarantee sporicidal sterilization.



The CD Check Strips can be placed throughout the area to be decontaminated to demonstrate that the chlorine dioxide gas reached all areas. The color-changed strips can be kept as part of your facility log to provide documented evidence that the decontamination process was performed.

The colored block changes color from a purple to a blue tone when exposed to chlorine dioxide gas.



To use, simply place the strips in the area to be decontaminated and then compare the color of the exposed strip to the color of a new strip to detect the color change. CD Check Strips come at a fraction of the cost of Biological Indicators (BI's), and without incubation and analysis of BI's, providing an easy method to document your gaseous chlorine dioxide decontamination process.

Product Number	Description
CHK-001	1 package/200 strips



50 Tannery Rd, Suite 1. Branchburg, NJ 08876 Tel:(908) 236-4100 Fax:(908) 236-2222

www.clordisys.com

31-030 Rev D